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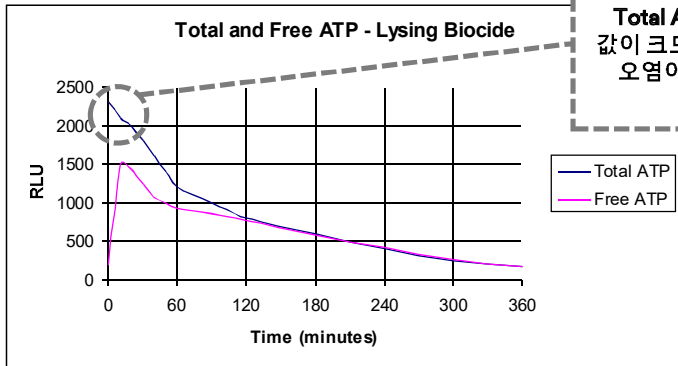
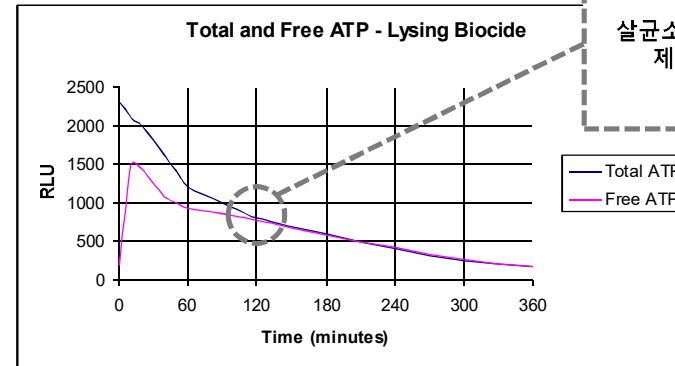
Case Study: Quat / Gluteraldehyde Treatment of a Dirty Water System

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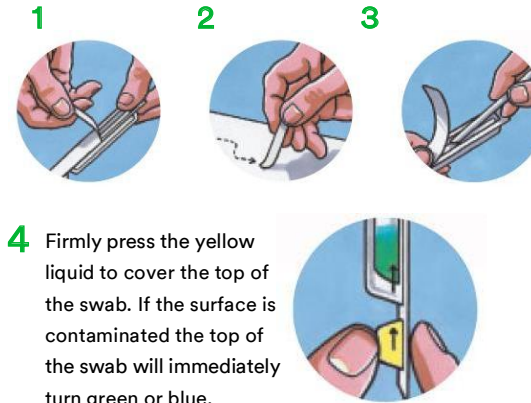
3M Health Care Academy

© 3M 2015. All Rights Reserved

3M Clean-Trace Surface Protein Instant

4 Firmly press the yellow liquid to cover the top of the swab. If the surface is contaminated the top of the swab will immediately turn green or blue.

3M Health Care Academy

© 3M 2015. All Rights Reserved

3M Protein Swabs

	3M Clean-Trace Surface Protein Allergen	3M Clean-Trace Surface Protein Plus	3M Clean-Trace Surface Protein Instant
Product Code	ALLTEC60	PRO50 and PRO100	CIT50
Detects	Protein, reducing sugars & agents	Protein, reducing sugars & agents	Protein only
Application	Validated for a range of allergenic proteins, including egg, milk, gluten, soy and peanut	General protein detection	Simpler & cheaper than PRO50. CIT50 is ideal for smaller food organizations where price is an issue.
Protocol	55°C for 15min	Room temperature for 10min	Room temperature for a few seconds
Limit of detection (using BSA)	3µg protein	50µg protein	50µg protein
Quantification	Semi-quantitative with 4 levels	Semi-quantitative with 4 levels	No, just Clean/Dirty
Handle colour	White	Purple	No handle
Swab tip	High-sensitivity swab tip	Regular foam tip	No swab tip

3M Health Care Academy

© 3M 2015. All Rights Reserved